


DinoDirect Installation Notice for Android Devices

The DinoDirect app is designed to connect Dino-Lite devices directly to Android devices via OTG. It was developed using an API version that ensures essential camera functionality and compatibility with Dino-Lite microscopes. However, due to changes in Google's policies regarding camera device usage in newer Android versions, the app cannot be listed on Google Play. To provide access to DinoDirect, the app is available as a downloadable APK.

Please follow the steps below to download and install the DinoDirect app:

1. Download the APK version of DinoDirect only from the official Dino-Lite website (www.dino-lite.com) for security reasons. If the installation does not execute automatically, please proceed with the steps below.
2. Tap "**More details**" on the pop-up message.

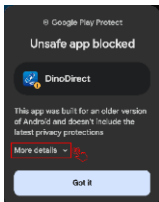

3. Select "**Install anyway**" to start the installation. The installation should then proceed automatically.

